

How to generate calibration and control reports from the Lx Maintenance software...

- 1) Goto **Start** → **Programs** → **Luminex** → **LXR** → **Lx Maintenance**
- 2) Under the **Tools** menu, select **Log Reports**
- 3) Under **Report Selection**, select **Calibration History**
- 4) Under **Date Range**, set the range of dates and times to include the reports desired
- 5) Leave the default setting in the **Success** and **Lot Numbers** fields unless you need a specific report file to be generated
- 6) Under **Phase**, choose the desired calibration/control beads for the report
- 7) Press the **View Button**
- 8) To save the report, click on the **File** menu and select **Save Report**
- 9) To print the report, click on the **File** menu and select **Print**
- 10) For **Control**, select **Control History** under **Report Selection** and repeat steps 4-9