

How do I generate the system log files from the Luminex Lx Maintenance software...

- 1) Goto **Start** → **Programs** → **Luminex** → **LXR** → **Lx Maintenance**
- 2) Under the **Tools** menu, select **Log Reports**
- 3) Under **Report Selection**, select **Log Messages**
- 4) Under **Date Range**, make sure you include the date of the last time it failed as well as a couple of days ahead and before the failure
- 5) Under **Message Type**, please select **All Message Types**
- 6) Press **View**
- 7) Under the **File** menu, select **Save Report**